



Instructions
Gebrauchsanleitung
Instrucciones
Istruizione

FLEXI **TESTER**
FTS300

Hydrogen Peroxide
(5 - 50 mg/l)




WATER TESTING
MADE IN GERMANY

Reagents
Reagenzien
Réactifs
Reactivos
Reagenti



FTS300

TbsHAFF
Acidifying PT Rapid
TbsRHP
Hyd. Peroxide Rapid

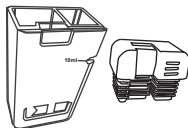

no detergent!



clean properly



WATER TESTING
MADE IN GERMANY

(GB)

CAUTION!

Reagents for water-analysis only!
Do not eat! Keep away from children!
Store cool and dry!
Clean tester properly before testing!

(F)

Attention!

Les pastilles sont exclusivement destinées aux analyses chimiques. NE PAS AVALER. Tenir hors de portée des enfants. Conserver dans un endroit frais et sec. Bien rincer le testeur avant chaque utilisation!

(E)

¡Atención!

Pastillas para el análisis de aguas, solamente para análisis químicos!
No para tomar!
No debe llegar a las manos de niños!
Consérvese en lugar fresco y seco!
Limpiar el aparato antes de usar!

(I)

Attenzione!

Reagenti unicamente per analisi acqua!
Non ingerire!
Tenere fuori dalla portata dei bambini!
Mantenere in luogo asciutto e fresco
Dopo l'utilizzo pulire ed asciugare accuratamente il tester.

(DE)


Achtung!

Reagenzien nur für chemische Analysen!
Nicht einnehmen! Darf nicht in die Hände von Kindern gelangen! Kühl und trocken lagern!




1

Fill with 10ml of
water to be tested

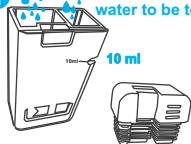


2

1 x Acidifying PT
Rapid
1 x +Hyd. Peroxide
HR Rapid

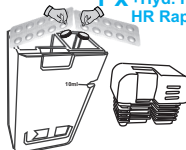

3

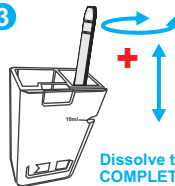

Dissolve tablet
COMPLETELY!

4

2 min.



5



6: Insert FTS300 stick
Hyd. Peroxide
5 - 50 mg/l

no detergent!



clean properly

