



- H315 Causes skin irritation.
- H319 Causes serious eye irritation.
- P101 If medical advice is needed, have product container or label at hand.
- P102 Keep out of reach of children.
- P264 Wash with water thoroughly after handling.
- P270 Do no eat, drink or smoke when using this product.
- P280 Wear protective gloves/protective clothing/eye protection/face protection.
- P302+352 IF ON SKIN: Wash with plenty of soap and water.
- P305+351+338 IF IN EYES: Rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing.
- P321 Specific treatment (see on this label).
- P332+313 If eye irritation persists: Get medical advice/attention.
- P337+313 If skin irritation occurs: Get medical advice/attention.
- P362 Take off contaminated clothing and wash before reuse.

NOTRUF/EMERGENCY: +49 (0)89-19240 (24h)



Total Hardness

Total Hardness

Tablet-count method

Gesamthärte

Tabletten-Zählverfahren

Dureté totale

Methode de comptage à pilule

Durezza totale

Metodo di conteggio de pastiglie

Dureza total

Método Contador de tabletas

Dureza total


Método de contagem de comprimidos

Total hårdhet

Tablettäggelings-metoden


Totale hardheid

färbet-telmethode

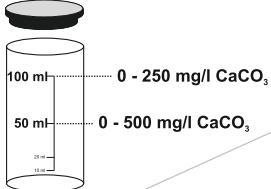


1


Total Hardness

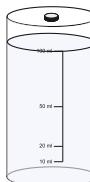

2

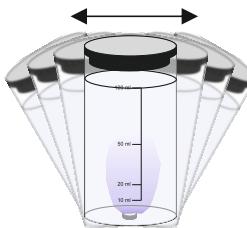


3

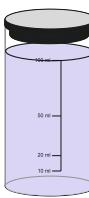

4




5

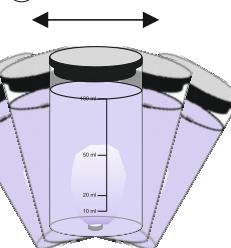4.5
4.5
4.5
4.5

Repeat until

=

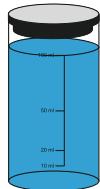

$$50 \text{ ml: } (\text{ } \times 40) - 20 = \text{mg/l CaCO}_3$$

$$100 \text{ ml: } (\text{ } \times 20) - 10 = \text{mg/l CaCO}_3$$



	mmol/l	mval/l	°dH	°e	°f
1 mg/l CaCO ₃	0.01	0.02	0.056	0.07	0.1

Alkaline Earth Ions
Erdalkali-Ionen
Ions alcalino-terreux
Ioni alcalino terrosi
Iones alcalina terrosos
Lões alcalino-terrosos
Ioner fra alkalsk jord
aardalkali-ionen